


---

# MicroPulser Electroporator

## Instruction Manual and Applications Guide

Catalog #1652100

For technical service call your local Bio-Rad office or in the U.S. call 1-800-4BIORAD (1-800-424-6723). On the Web at [bio-rad.com](http://bio-rad.com).

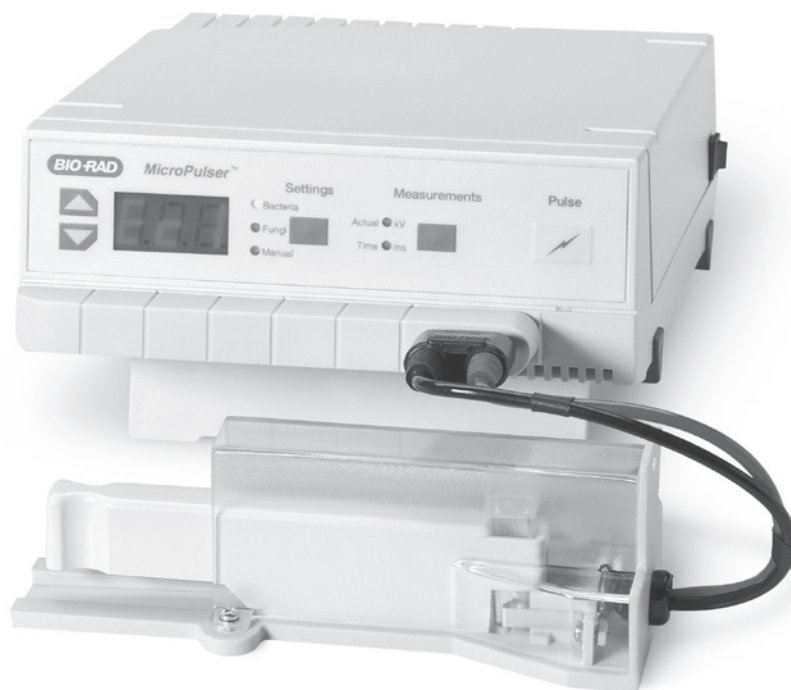

**BIO-RAD**

# Warranty

Bio-Rad Laboratories warrants the MicroPulser Electroporator against defects in materials and workmanship for 1 year. If any defects occur in the instrument during this warranty period, Bio-Rad Laboratories, Inc. will, at Bio-Rad's option, repair or replace the defective parts free of charge. The following defects, however, are specifically excluded:

1. Defects caused by improper operation.
2. Repair or modification done by anyone other than Bio-Rad Laboratories or an authorized agent.
3. Use of fittings or other spare parts supplied by anyone other than Bio-Rad Laboratories.
4. Damage caused by accident or misuse.
5. Damage caused by disaster.
6. Corrosion due to use of improper solvent or sample.

For any inquiry or request for repair service, contact Bio-Rad Laboratories after confirming the model, serial number, invoice number, and purchase order number of your instrument.

Model: Micropulser Electroporator \_\_\_\_\_

Catalog No.: \_\_\_\_\_

Date of Delivery: \_\_\_\_\_

Serial No.: \_\_\_\_\_

Invoice No.: \_\_\_\_\_

Purchase Order No.: \_\_\_\_\_

# Table of Contents

<b>Section 1</b>	<b>Safety Information</b> . . . . .	<b>1</b>
	1.1 Electrical Hazards . . . . .	1
	1.2 Mechanical Hazards . . . . .	1
	1.3 Other Safety Precautions . . . . .	1
<b>Section 2</b>	<b>Introduction</b> . . . . .	<b>2</b>
	2.1 Overview of Electroporation Theory . . . . .	2
	2.2 Manipulation of Instrument Parameters . . . . .	4
<b>Section 3</b>	<b>Factors Affecting Electroporation</b> . . . . .	<b>5</b>
	3.1 Cell Growth . . . . .	5
	3.2 DNA . . . . .	5
	3.3 Electroporation Media . . . . .	6
<b>Section 4</b>	<b>MicroPulser Electroporator Operating Instructions</b> . . . . .	<b>8</b>
	4.1 Setting up the MicroPulser Electroporator . . . . .	8
	4.2 Operation of the MicroPulser Electroporator . . . . .	8
	4.2.1 Selecting Preprogrammed Settings . . . . .	8
	4.2.2 Using the MicroPulser Electroporator in Manual Mode . . . . .	9
	4.2.2A Changing the Voltage . . . . .	9
	4.2.2B Truncating the Pulse . . . . .	9
	4.2.3 Pulse Function . . . . .	10
	4.2.4 Measurements . . . . .	10
	4.3 Electroporation Using the MicroPulser Electroporator . . . . .	10
<b>Section 5</b>	<b>High-Efficiency Electrotransformation of <i>Escherichia coli</i></b> . . . . .	<b>11</b>
	5.1 Preparation of Electrocompetent Cells . . . . .	11
	5.2 Electroporation . . . . .	11
	5.3 Solutions and Reagents for Electroporation . . . . .	12

<b>Section 6</b>	<b>Electroporation of <i>Staphylococcus aureus</i></b> .....	<b>12</b>
	6.1 Preparation of Electrocompetent Cells. ....	12
	6.2 Electroporation .....	13
	6.3 Solutions and Reagents for Electroporation .....	13
<b>Section 7</b>	<b>Electroporation of <i>Agrobacterium tumefaciens</i></b> .....	<b>14</b>
	7.1 Preparation of Electrocompetent Cells. ....	14
	7.2 Electroporation .....	14
	7.3 Solutions and Reagents for Electroporation. ....	15
<b>Section 8</b>	<b>Electroporation of <i>Saccharomyces cerevisiae</i></b> .....	<b>15</b>
	8.1 Preparation of Electrocompetent Cells. ....	15
	8.2 Electroporation .....	16
	8.3 Solutions and Reagents for Electroporation .....	16
<b>Section 9</b>	<b>Electroporation of <i>Schizosaccharomyces pombe</i></b> .....	<b>16</b>
	9.1 Preparation of Electrocompetent Cells. ....	16
	9.2 Electroporation .....	17
	9.3 Solutions and Reagents for Electroporation. ....	17
<b>Section 10</b>	<b>Electroporation of <i>Dictyostelium discoideum</i></b> .....	<b>18</b>
	10.1 Preparation of Electrocompetent Cells. ....	18
	10.2 Electroporation .....	18
	10.3 Solutions and Reagents for Electroporation .....	19
<b>Section 11</b>	<b>Electroporation of <i>Pichia pastoris</i></b> .....	<b>19</b>
	11.1 Preparation of Electrocompetent Cells .....	19
	11.2 Electroporation. ....	20
	11.3 Solutions and Reagents for Electroporation .....	20
<b>Appendix I</b>	<b>References</b> .....	<b>21</b>
<b>Appendix II</b>	<b>Troubleshooting Guide for the MicroPulser Electroporator</b> .....	<b>22</b>
<b>Appendix III</b>	<b>Product Information</b> .....	<b>25</b>

# Section 1

## Safety Information

### Read This Information Carefully Before Using the MicroPulser Electroporator.

The MicroPulser Electroporator meets the safety requirements of EN 61010-1 and the electromagnetic compatibility (EMC) requirements of EN 61326 (for Class B, including flicker and harmonics).

### 1.1 Electrical Hazards

The MicroPulser Electroporator produces voltages up to 3,000 V and can pass very high currents. When charged to maximum voltage, the instrument stores about 50 J. A certain degree of respect is required for energy levels of this order. Safety system features prevent operator access to the recessed input jacks and to the recessed electrode contacts inside the sample chamber. These mechanical interlocks should never be circumvented.

There is high voltage present whenever the yellow pulse button is depressed and “PLS” is shown in the light emitting diode display on the front of the instrument. If the capacitor has been partially charged but not fired (for example, when the charging cycle has been interrupted before the pulse is delivered), some charge may remain on the internal capacitor. However, the user cannot make contact due to the system safety features.

### 1.2 Mechanical Hazards

The MicroPulser Electroporator contains a patented arc-protection circuit that dramatically reduces the incidence of arcing in the cuvette when high voltage is delivered into the sample. The unit incorporates a circuit which senses the beginning of an arc and diverts current from the sample within ~5  $\mu$ s, preventing or greatly reducing mechanical, visual, and auditory phenomena at the shocking chamber. Should an arc occur, the sample chamber is effective in containing these small discharges; nonetheless, we strongly recommend wearing safety glasses when using the instrument.

Do not use the MicroPulser Electroporator with samples suspended in conductive media (refer to Section 3.3 for information on sample resistance).

### 1.3 Other Safety Precautions

- Turn the unit off when not attended
- Avoid spilling any liquids onto the apparatus. Use only a paper towel or a cloth wet with either water or alcohol to clean the outside surfaces of the MicroPulser Electroporator
- Use only the Bio-Rad cables supplied with the MicroPulser Electroporator
- Only use the shocking chamber in the assembled condition. Do not attempt to circumvent the protection of the shocking chamber or use it while disassembled
- Verify the display segments periodically
- Do not use the MicroPulser Electroporator if obvious case damage exists that exposes part of the inside of the unit
- No user-serviceable parts are contained within the MicroPulser Electroporator; the case should only be opened by properly trained personnel

**Warning:** The MicroPulser Electroporator generates, uses, and radiates radio frequency energy. If it is not used in accordance with the instructions given in this manual, it may cause interference with radio communications. The MicroPulser Electroporator has been tested and found to comply with the limits for Class A computing devices (pursuant to Subpart J of Part 15 of FCC Rules), which provide reasonable protection against such interference when operated in a commercial environment. Operation of this equipment in a residential area is likely to cause interference. In this case the user will be required, at their own expense, to take whatever measure may be required to correct the interference.

## Section 2 Introduction

### 2.1 Overview of Electroporation Theory

The MicroPulser Electroporator is used for the electroporation of bacteria, yeast, and other microorganisms, where a high-voltage electrical pulse is applied to a sample suspended in a small volume of high-resistance medium. The system consists of a pulse generator module, a shocking chamber, and a cuvette with incorporated electrodes (Figure 1). The sample is placed between the electrodes in the cuvette. The MicroPulser Electroporator module contains a capacitor, which is charged to a high voltage; the module then discharges the current in the capacitor into the sample in the cuvette.

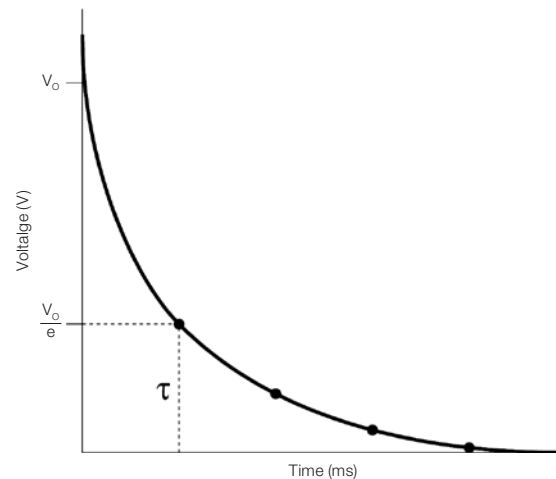


Fig. 1. MicroPulser Electroporator consists of pulse generator module, shocking chamber, and cuvette.

The capacitance discharge circuit of the MicroPulser Electroporator generates an electrical pulse with an exponential decay waveform (Figure 2). When the capacitor is discharged into the sample, the voltage across the electrodes rises rapidly to the peak voltage (also known as the initial voltage,  $V_0$ ) and declines over time,  $t$ , as follows:

$$V_t = V_0 [e^{-(t/\tau)}],$$

where  $\tau = R \times C$ , the time constant, a convenient expression of the pulse length. The resistance of the circuit,  $R$ , is expressed in ohms, and the capacitance of the apparatus,  $C$ , is expressed in microfarads. According to the equation above,  $\tau$  is the time over which the voltage declines to  $1/e$  (~37%) of the peak value. The internal circuitry of the MicroPulser Electroporator is designed to provide optimum electroporation of *Escherichia coli* and *Saccharomyces cerevisiae*, as well as many other microorganisms, in which the optimum transformation efficiency occurs at a time constant of approximately 5 ms. These electroporation conditions are achieved by using a 10 mF capacitor and by placing a 600  $\Omega$  resistor in parallel with the sample cuvette along with a 30  $\Omega$  resistor in series with the sample cuvette.



**Fig. 2. Exponential decay pulse from a capacitance discharge system.** When the capacitor, charged to an initial voltage,  $V_0$ , is discharged into cells, the voltage applied to the cells decreases over time so that at time  $t = \tau$ , the voltage is  $(1/e) \times V_0$  of the initial value.

In addition to the time constant, the electric field strength is the other instrument parameter that is important in determining transformation efficiency. The electric field strength,  $E$ , is the voltage applied between the electrodes and is described by

$$E = V / d,$$

where  $V$  is the voltage applied and  $d$  is the distance (cm) between the electrodes. The strength of the electric field and the size of the cells determine the voltage drop across each cell, and it is this voltage drop that may be the important manifestation of the voltage effect in electroporation.

The purpose of the 30  $\Omega$  series resistor in the MicroPulser Electroporator is to protect the instrument circuitry should arcing occur. Under normal operation, when samples are in high-resistance media, this resistor will not affect the voltage applied to the sample. However, this resistor will significantly decrease the voltage applied to the sample if the resistance of the sample is low. The fractional drop in voltage applied to the sample is given by

$$R_{30} / (R_{30} + R_{\text{sample}}).$$

When  $R_{\text{sample}}$  is 600  $\Omega$ , there is a 5% voltage drop to the sample:  $30 / (30 + 600) = 0.048$ .

For this reason, electroporation with the MicroPulser Electroporator should not be performed in solutions with a resistance of less than ~600  $\Omega$ , including samples in which the growth medium was not adequately removed from the cells, DNA samples containing salt contributed by residual sodium chloride, or ligation mixtures. The MicroPulser Electroporator can measure the resistance of the sample and will not pulse into very low-resistance media.


## 2.2 Manipulation of Instrument Parameters

Several parameters on the MicroPulser Electroporator may be altered to achieve maximum transformation efficiency. These include the field strength,  $E$ , the time constant,  $\tau$ , and the width of a truncated exponential decay pulse. The field strength may be manipulated in two ways. First, voltages between 200 and 3,000 V may be set directly on the MicroPulser Electroporator. This parameter is the most easily controlled.

The process of varying the voltage while keeping all other conditions unchanged is the basis for most electroporation optimization procedures. Second, using cuvettes with different electrode gap widths permits a means of changing the field strength. For electroporation of microorganisms, 0.1 and 0.2 cm gap cuvettes are most often used. Electroporation of *E. coli* is generally carried out at a voltage of 1.8 kV ( $E = 18$  kV/cm) when electroporating cells in 0.1 cm cuvettes and at a voltage of 2.5 kV ( $E = 12.5$  kV/cm) when electroporating cells in 0.2 cm cuvettes. These electroporation conditions are preprogrammed into the MicroPulser Electroporator as programs Ec1 ( $V = 1.8$  kV) and Ec2 ( $V = 2.5$  kV) in the bacterial settings menu. In addition, a third program, Ec3 in the bacterial settings menu, delivers a voltage of 3.0 kV ( $E = 15$  kV/cm in 0.2 cm cuvettes), which we have found results in even higher transformation efficiency compared to electroporation at 2.5 kV.

The time constant may be altered by changing the sample resistance. The sample resistance may be manipulated in two ways. First, increasing the salt or buffer concentration of the electroporation medium decreases the resistance of the sample and vice versa, resulting in a change in the time constant. Second, the volume of the sample in the cuvette is inversely proportional to the resistance of the sample; decreasing the sample volume increases the sample resistance. This effect of volume on sample resistance is most noticeable in low-resistance media. These effects are discussed further in Section 3.3.

The MicroPulser Electroporator also includes a means to truncate the exponential decay pulse sooner than the expected time constant as long as the voltage is greater than 600 V. When the pulse is terminated by the MicroPulser Electroporator, voltage is applied to the sample only for the specified time, which may be between 1.0 and 4.0 ms. Figure 3 shows how this waveform differs from the true exponential decay pulse.

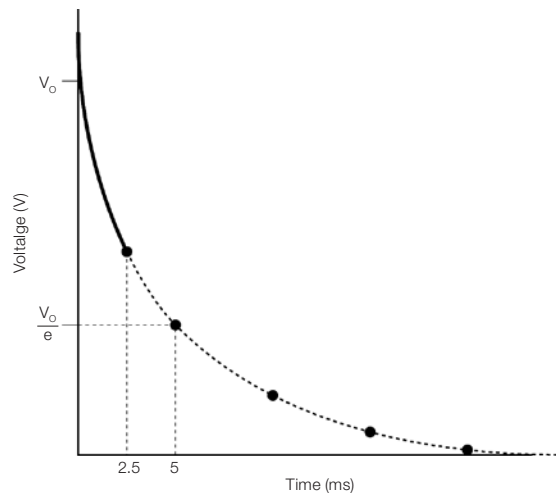

**Fig. 3. Truncation of an exponential decay pulse by the MicroPulser Electroporator.** The solid line shows the voltage applied to the cells as a function of time during a pulse terminated after 2.5 ms. The dashed line shows the voltage that would normally be applied to the cells during a true exponential decay pulse.



## Section 3

# Factors Affecting Electroporation

The electrical conditions for the electroporation of microorganisms have been verified through years of research (see Chang et al. 1992 and Nickoloff 1995 for overviews as well as for protocols on electroporation of numerous species). For many microorganisms, optimum electrotransformation occurs under electrical conditions relatively similar to those used for *E. coli* and *S. cerevisiae*, two species that are most commonly used in research today. For electroporation of *E. coli*, conditions reported as being used most often are 0.2 cm cuvettes containing 40  $\mu$ l of cells at a voltage of 2.5 kV and a time constant of  $\sim$ 5 ms. For electroporation of *S. cerevisiae*, conditions reported as being used most often are 0.2 cm cuvettes containing 40  $\mu$ l of cells at a voltage of 1.5 kV and a time constant of  $\sim$ 5 ms. For many bacteria, including species of *Salmonella*, *Pseudomonas*, *Helicobacter*, *Borrelia*, *Streptococcus*, *Lactococcus*, and *Enterococcus*, the conditions for electroporation are identical to those used for *E. coli*. For many other bacterial species, altering the field strength will often result in higher electrotransformation efficiency. A similar case is found with other species of yeast.

The MicroPulser Electroporator is designed to precisely deliver those pulse parameters needed for the highest transformation efficiency of *E. coli* and *S. cerevisiae*. The time constant has been set at 5 milliseconds when working with high-resistance samples. For these organisms, the MicroPulser Electroporator has preprogrammed settings for delivery of the correct voltage when electroporating *E. coli* in either 0.1 or 0.2 cm cuvettes, or when electroporating *S. cerevisiae* in either 0.2 or 0.4 cm cuvettes.

### 3.1 Cell Growth

For most bacterial species, the highest transformation efficiencies are obtained when cells are harvested in early- to mid-log growth. For *E. coli*, as the cells reach stationary phase, the transformation efficiency will decline precipitously (Dower 1990). In contrast, most yeast species are generally harvested in mid- to late-log growth. For *S. cerevisiae*, the transformation efficiency increases as much as 60-fold from early- to late-log cultures (Becker and Guarente 1991). The optimal portion of the growth phase to harvest cells is generally dependent on the cell type. When preparing competent cells of a new species, it is generally best to employ conditions worked out for use with the same genus. Suggestions for factors to consider and general methods for producing electrocompetent cells are discussed in the articles by Dower et al. 1992 and Trevors et al. 1992.

### 3.2 DNA

While the majority of electroporation applications involve delivery of plasmid DNA to cells, it should be mentioned that nearly any type of molecule can be introduced into cells by electroporation, including RNA, proteins, carbohydrates, and small molecules. With few exceptions, when delivering autonomously replicating plasmids, the highest transformation efficiencies are obtained when electroporating supercoiled plasmid. However, electroporating plasmid that will integrate into the host genome is usually most efficient using linear plasmid. For example, *Candida*, *Pichia*, and *Tetrahymena* are transformed more efficiently when transformed with linearized rather than with supercoiled integrating plasmids.

In both *E. coli* and *Listeria monocytogenes*, the transformation efficiency of relaxed circular plasmid is only slightly lower than that of supercoiled plasmid (Leonardo and Sedivy 1990; Park and Stewart 1990). However, linear plasmid is about  $10^3$ – $10^4$ -fold less efficient than the corresponding circular plasmid in both *E. coli* and *Streptococcus pyogenes* (Shigekawa and Dower 1988; Simon and Ferretti 1991). Electroporation efficiency per mole of plasmid generally decreases as the plasmid size increases in numerous species, including *E. coli* (Leonardo and Sedivy 1990; Siguret et al. 1994), *Pseudomonas aeruginosa* (Dennis and Sokol 1995), and *Streptococcus thermophilus* (Somkuti and Steinberg 1988). However, in some species, including *Lactococcus lactis* (Holo and Nes 1995), *Enterococcus faecalis* (Cruz-Rodz and Gilmore 1990), and *Clostridium perfringens* (Allen and Blaschek 1990), transformation efficiency appears to be independent of plasmid size up to 20–30 kb.


Although transformation of most microorganisms has been accomplished using plasmid DNA isolated by a variety of methods, the plasmid purity has an effect on transformation efficiency. Significantly lower transformation efficiencies are generated with unpurified miniprep plasmid DNA than with plasmid DNA purified by a variety of procedures. Plasmid produced using the Bio-Rad Quantum matrix is as efficient as CsCl-purified plasmid for transformation of microorganisms.

Generally, for all types of microorganisms, the frequency of transformation increases with increasing DNA concentration in the electroporation buffer. For *E. coli*, the frequency of transformation (transformants/survivor) is dependent on DNA concentration over at least six orders of magnitude (10 pg/ml to 7.5  $\mu$ g/ml); within this range the DNA concentration determines the probability that a cell will be transformed. At the higher DNA concentrations, up to 80% of the survivors are transformed (Dower et al. 1988). Because the number of transformants recovered is the product of the transformation frequency and the number of cells present, the transformation efficiency (transformants/ $\mu$ g DNA) increases with cell concentration over the range of  $10^9$  to at least  $3 \times 10^{10}$  cells/ml. Therefore, to obtain a high transformation frequency, use high DNA concentration. To obtain high transformation efficiency, use high cell concentration and low DNA concentration to avoid cotransformations. In each case, a small sample volume (20–50  $\mu$ l) allows economical use of DNA and cells (see Dower et al. 1988 for a detailed discussion of these factors).

### 3.3 Electroporation Media

The MicroPulser Electroporator is designed for use with samples in high-resistance media (>600  $\Omega$ ). For this reason, when preparing electrocompetent cells, it is important to wash cells thoroughly to remove all traces of growth medium. Failure to thoroughly remove the growth medium from the cells may result in the sample arcing during electroporation. Cells should be washed at least three times with water or with nonionic solutions, such as glucose, glycerol, sucrose, sorbitol, or polyethylene glycol. For many microorganisms, glycerol is a convenient electroporation medium, since it is recommended as a cryoprotectant for storage of cell cultures.

Figures 4A and B show the effect of concentration of several biologically important ionic solutions on sample resistance. **Note:** (1) volume has a significant effect on sample resistance for ionic solutions, and sample resistance is inversely proportional to the volume of solution in the cuvette; (2) the resistance of a solution containing divalent ions is lower than that of a solution containing the same concentration of monovalent ions; (3) the resistance of a buffered solution is affected by its pH.

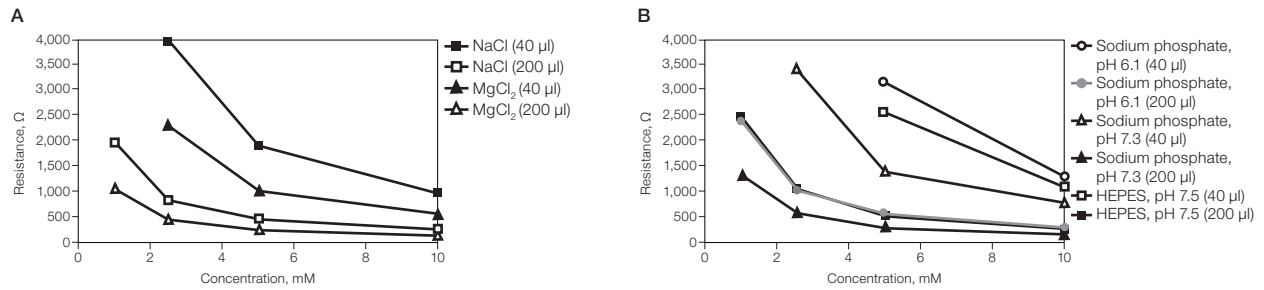

**Fig. 4. Resistance of A, solutions of NaCl and MgCl<sub>2</sub> and B, buffers of sodium phosphate at pH 6.1 and 7.3 and HEPES at pH 7.5.** Resistance was measured in 0.2 cm cuvettes containing either 40 or 200 µl of solution at room temperature.

Even small concentrations of ionic compounds significantly reduce the resistance of the sample and may cause arcing. Residual salt from ethanol precipitation of DNA should be reduced by washing the DNA pellet prior to resuspending it in either water or Tris-EDTA (TE). Table 1 shows that although adding a solution of plasmid in 10 mM Tris, pH 8.0, 1 mM EDTA to water does reduce the sample resistance, this should not result in the inability to electroporate a sample in the MicroPulser Electroporator. DNA may be used directly from enzyme reactions for transformation, but the final salt concentration in the electroporation sample should be kept below ~5 meq for high-voltage operation. Finally, ligation mixtures may be used for transformation, but only in very low quantities or when the ionic strength is reduced by dilution (Willson and Gough 1988), dialysis (Heery and Dunican 1989; Jacobs et al. 1990), or ethanol precipitation (Böttger 1988; Zabarovsky and Winberg 1990).

**Table 1. Resistance of water in 0.2 cm cuvettes to which TE has been added.\***

Sample	R <sub>sample</sub> * Ω	
	(40 µl volume)	(200 µl volume)
Water	>6 x 10 <sup>5</sup>	>6 x 10 <sup>5</sup>
Water + 1 µl TE	>6 x 10 <sup>5</sup>	35,000
Water + 5 µl TE	11,200	8,700
Water + 10 µl TE	4,850	4,700

\* The resistance of 0.2 cm cuvettes containing either 40 or 200 µl water and the indicated volume of TE (10 mM Tris, pH 8.0, 1 mM EDTA) was measured at 1,000 V.

## Section 4

# MicroPulser Electroporator Operating Instructions

Refer to Figure 1 for a view of the components of the MicroPulser Electroporator and to Figure 5 for a schematic of the buttons and light emitting diodes (LEDs).

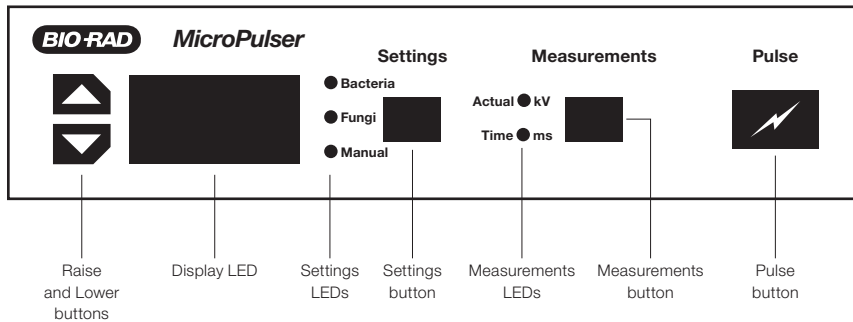


Fig. 5. MicroPulser Electroporator control panel.

### 4.1 Setting up the MicroPulser Electroporator

1. Connect the black power cord to the rear panel of the MicroPulser pulse generator module. Plug the cord into a wall outlet or power strip.
2. Pull down the fold-down foot on the underside of the MicroPulser Electroporator. Insert this foot into the track on the base of the shocking chamber. Insert the shocking chamber slide into the shocking chamber.
3. Connect the leads from the shocking chamber to the output jacks on the front panel of the MicroPulser Electroporator; polarity is not important to the electroporation process.
4. Turn on the apparatus using the power switch on the right rear panel. The LED display should illuminate and read "Ec1," and the LED next to the Bacteria Settings should be illuminated.

### 4.2 Operation of the MicroPulser Electroporator

#### 4.2.1 Selecting Preprogrammed Settings

The MicroPulser Electroporator is preprogrammed with settings for electroporation of several commonly used organisms.

Bacteria Settings program:

Mnemonic	Organism	Cuvette size, cm	Parameters		
			Voltage, kV	Number of pulses	Time constant, ms
Ec1	<i>E. coli</i>	0.1	1.8	1	—
Ec2	<i>E. coli</i>	0.2	2.5	1	—
StA	<i>S. aureus</i>	0.2	1.8	1	2.5
Agr	<i>A. tumefaciens</i>	0.1	2.2	1	—
Ec3	<i>E. coli</i>	0.2	3.0	1	—

Fungi Settings program:

Mnemonic	Organism	Cuvette size, cm	Parameters		
			Voltage, kV	Number of pulses	Time constant, ms
Sc2	<i>S. cerevisiae</i>	0.2	1.5	1	—
Sc4	<i>S. cerevisiae</i>	0.2	3.0	1	—
ShS	<i>S. pombe</i>	0.2	2.0	1	—
Dic	<i>D. discoideum</i>	0.4	1.0	2	1.0
Pic	<i>P. pastoris</i>	0.2	2.0	1	—

Pressing the “Settings” button cycles the Settings LED between “Bacteria,” “Fungi,” and “Manual.” To change from one bacterium setting to another, while the Settings LED next to Bacteria is lit, press the “Raise” and “Lower” buttons to the left of the display LED to cycle between the different programs. When the mnemonic is displayed, the parameters associated with the mnemonic are automatically selected.

When the LED next to Fungi is lit, the mnemonic for the fungi programs is displayed. Press the “Raise” and “Lower” buttons to the left of the display to cycle between the different programs. When the mnemonic is displayed, the parameters associated with the mnemonic are automatically selected.

While a program mnemonic is displayed for either a Bacteria or Fungi Setting, simultaneously pressing both the “Raise” and “Lower” buttons shows the program parameters selected on the display LED. The display LED first shows the voltage value, then displays a “t” followed by the time in milliseconds, then, if time and multiple pulses are associated with a program, displays “P” followed by “2,” indicating that two successive pulses are given. If no “t” is given, the pulse is not truncated, and if no “P” is given, there is a single pulse.

## 4.2.2 Using the MicroPulser Electroporator in Manual Mode

### 4.2.2A Changing the Voltage

Press the “Settings” button to illuminate the LED next to “Manual.” The display LED now shows the voltage (in kV). Pressing the “Raise” and “Lower” buttons to the left of the display LED allows selection of the desired voltage in the range from 0.20 to 3.00 kV. If the instrument was just turned on, the display LED shows “0.00.”

### 4.2.2B Truncating the Pulse

Press both “Raise” and “Lower” buttons simultaneously while the “Manual” LED is illuminated. The display LED now shows “t—” and indicates the pulse time that has been selected for the pulse. The default setting when the power is turned on is the standard exponential decay pulse, or no pulse truncation, which is indicated by the two dashes. Releasing only the “Lower” button results in the display LED showing the time of the truncated pulse in milliseconds. The display LED initially changes to “t1.0” and rises in 0.1 ms increments to “t4.0.” This permits truncating the exponential decay pulse between 1 and 4 ms. Simultaneously pressing both “Raise” and “Lower” buttons, then releasing only the “Raise” button, results in lowering the indicated truncation time.

### 4.2.3 Pulse Function

Pressing the “Pulse” button results in the capacitor charging to the set voltage; during this time “PLS” is shown on the display LED. A tone will sound to indicate that the pulse has been delivered. When multiple pulses are delivered by one of the built-in programs, “PLS” is shown on the display LED during the entire time and a tone sounds each time a pulse is delivered. To manually deliver multiple pulses, after the tone sounds from the first pulse, press the pulse button again.

If a lower pitched tone sounds, accompanied by “Arc” being shown on the display LED, the arc prevention and quenching (ARQ) system has been actuated and the pulse has been terminated. This is usually an indication of attempted cuvette arcover, but may also occur if the sample resistance is too low. Since the energy delivered during such an ARQ event is low, it is usually possible to pulse the sample again at parameters which will not result in an arc and still produce acceptable results. However, it is not advisable to use a sample in which two arc events have occurred.

### 4.2.4 Measurements

Pressing the “Measurements” button results in illumination of the “Actual kV” LED. This indicates that the display LED shows the actual voltage delivered (in kV) during the last pulse. If the instrument was just turned on and no pulse has been given, the display LED shows “0.00.” Pressing the “Measurements” button again results in illumination of the “Time ms” LED. This indicates that the display LED shows the time constant (in ms) of the last pulse. If the instrument was just turned on and no pulse has been given, the display LED shows “0.00.” Holding the “Measurements” button toggles the display LED between the Actual voltage and the Time constant.

## 4.3 Electroporation Using the MicroPulser Electroporator

1. Place the cell suspension in an electroporation cuvette and tap the liquid to the bottom of the cuvette. Up to 0.4 ml (400  $\mu$ l) of solution may be placed in the 0.2 cm cuvette, and up to 80  $\mu$ l may be placed in the 0.1 cm cuvette. **Note:** Temperature may have a significant influence on transformation frequency. Electroporation of some organisms, including *E. coli* and *S. cerevisiae*, is more efficient in chilled cuvettes.
2. Insert the cuvette into the slide of the shocking chamber. Push the slide into the chamber until the cuvette makes firm contact with the chamber electrodes.
3. To charge the capacitor and deliver a pulse, press the yellow “Pulse” button; the display LED will show “PLS” until a tone sounds, indicating that the pulse has been given. The display LED will then show the program, the time constant, or the actual volts delivered, depending on the LED selected.
4. Withdraw the slide from the chamber, remove the cuvette, and process the sample.
5. The time constant and the actual voltage delivered to the sample are shown on the display LED by pressing the “Measurements” button. When the LED next to “Actual kV” is illuminated, the voltage is displayed in kilovolts. The time constant can be displayed by pressing the “Measurements” button again. The LED next to “Time ms” will be illuminated; the time constant is displayed in milliseconds.
6. To turn the unit off, press the power switch on the right rear panel. The sample chamber may now be safely disconnected, if desired. Never remove the sample chamber cover until the leads are disconnected.

## Section 5

# High-Efficiency Electrotransformation of *Escherichia coli*

Electroporation provides a method of transforming *E. coli* at efficiencies as high as  $10^9$  to  $10^{10}$  transformants/ $\mu\text{g}$ , which is greater than is possible with the best chemical methods. The following protocol describes a method for preparing and electrotransforming *E. coli* to high efficiencies.

### 5.1 Preparation of Electrocompetent Cells

See Ausubel et al. 1987 and Miller and Nickoloff 1995 for additional information.

1. Inoculate 500 ml of LB broth with 1/100 volume of a fresh overnight *E. coli* culture.
2. Grow the cells at 37°C shaking at 300 rpm to an OD 600 of ~0.5–0.7 (the best results are obtained with cells that are harvested at early- to mid-log phase; the appropriate cell density, therefore, depends on the strain and growth conditions).
3. Chill cells on ice for ~20 min. For all subsequent steps, keep the cells as close to 0°C as possible (in an ice-water bath) and chill all containers in ice before adding cells. To harvest, transfer the cells to a cold centrifuge bottle and spin at 4,000 x g for 15 min at 4°C.
4. Carefully pour off and discard the supernatant. It is better to sacrifice the yield by pouring off a few cells than to leave any supernatant behind.
5. Gently resuspend the pellet in 500 ml of ice-cold 10% glycerol. Centrifuge at 4,000 x g for 15 min at 4°C; carefully pour off and discard the supernatant.
6. Resuspend the pellet in 250 ml of ice-cold 10% glycerol. Centrifuge at 4,000 x g for 15 min at 4°C; carefully pour off and discard the supernatant.
7. Resuspend the pellet in ~20 ml of ice-cold 10% glycerol. Transfer to a 30 ml sterile Oakridge tube. Centrifuge at 4,000 x g for 15 min at 4°C; carefully pour off and discard the supernatant.
8. Resuspend the cell pellet in a final volume of 1–2 ml of ice-cold 10% glycerol. The cell concentration should be about  $1\text{--}3 \times 10^{10}$  cells/ml.

This suspension may be frozen in aliquots on dry ice and stored at  $-70^\circ\text{C}$ . The cells are stable for at least six months under these conditions.

### 5.2 Electroporation

1. Thaw the cells on ice. For each sample to be electroporated, place a 1.5 ml microfuge tube and either a 0.1 or 0.2 cm electroporation cuvette on ice.
2. In a cold, 1.5 ml polypropylene microfuge tube, mix 40  $\mu\text{l}$  of the cell suspension with 1 to 2  $\mu\text{l}$  of DNA (DNA should be in a low ionic strength buffer such as TE). Mix well and incubate on ice for ~1 min. (**Note:** It is best to mix the plasmids and cells in a microfuge tube since the narrow gap of the cuvettes prevents uniform mixing.)

3. Set the MicroPulser Electroporator to “Ec1” when using the 0.1 cm cuvettes. Set it to “Ec2” or “Ec3” when using the 0.2 cm cuvettes. See Section 4 for operating instructions.
4. Transfer the mixture of cells and DNA to a cold electroporation cuvette and tap the suspension to the bottom. Place the cuvette in the chamber slide. Push the slide into the chamber until the cuvette is seated between the contacts in the base of the chamber. Pulse once.
5. Remove the cuvette from the chamber and immediately add 1 ml of SOC medium to the cuvette. Quickly but gently resuspend the cells with a Pasteur pipet. The period between applying the pulse and transferring the cells to outgrowth medium is crucial for recovering *E. coli* transformants (Dower et al. 1988). Delaying this transfer by even 1 minute causes a threefold drop in transformation efficiency. This decline continues to a 20-fold drop by 10 minutes.
6. Transfer the cell suspension to a 17 x 100 mm polypropylene tube and incubate at 37°C for 1 hr, shaking at 225 rpm.
7. Check and record the pulse parameters. The time constant should be close to 5 milliseconds. The field strength can be calculated as actual kilovolts (kV) / cuvette gap (cm).
8. Plate on selective medium.

### 5.3 Solutions and Reagents for Electroporation

1. LB Broth: 10 g Bacto tryptone, 5 g Bacto yeast extract, 5 g NaCl; dissolve in 1.0 L water. Autoclave.
2. 10% (v/v) glycerol: 12.6 g glycerol (density = 1.26 g/ml) in 90 ml of water. Autoclave or filter sterilize.
3. TE: 10 mM Tris-HCl, pH 8.0, 1 mM EDTA.
4. SOC: 2% Bacto tryptone, 0.5% Bacto yeast extract, 10 mM NaCl, 2.5 mM KCl, 10 mM MgCl<sub>2</sub>, 10 mM MgSO<sub>4</sub>, 20 mM glucose.

## Section 6

# Electroporation of *Staphylococcus aureus*

### 6.1 Preparation of Electrocompetent Cells

See Lee 1995 for additional information.

1. Inoculate 3 ml of B2 broth in a 17 x 100 mm tube with a colony from a fresh *S. aureus* plate.
2. Incubate at 37°C overnight, shaking at 250 rpm.
3. Inoculate 1.5 ml of the overnight culture into 150 ml of fresh B2 broth in a 1 L flask. Incubate at 37°C, shaking at 250 rpm, to ~2 x 10<sup>8</sup> cells/ml. The doubling time of *S. aureus* is ~30 min at 37°C.
4. Chill the cells in an ice-water bath for 15 min to stop growth. Decant the cells into a sterile 500 ml centrifuge bottle. Harvest the cells by centrifugation at 12,000 x g for 15 min at 4°C.
5. Carefully pipet off the supernatant, keeping the cell pellet on ice.



6. Resuspend the cell pellet in 500 ml of sterile, ice-cold water. Pellet the cells by centrifugation at 12,000 x g for 15 min at 4°C; carefully remove the supernatant. Wash the cells two more times in 500 ml of sterile, ice-cold water.
7. Resuspend the cell pellet in 25 ml of sterile, ice-cold 10% glycerol. Transfer to a 30 ml sterile Oakridge tube. Pellet the cells by centrifugation at 12,000 x g for 15 min at 4°C; carefully remove the supernatant.
8. Resuspend the cell pellet in 2 ml of 10% glycerol; the final cell concentration should be  $\sim 1 \times 10^{10}$  cells/ml.
9. Dispense 250  $\mu$ l aliquots of the electrocompetent cells into sterile 1.5 ml microfuge tubes; freeze the cells in an isopropanol-dry ice bath, then store at  $-70^{\circ}\text{C}$ . The cells are stable for several months under these conditions.

## 6.2 Electroporation

1. Pipet the DNA samples (5 ng–2  $\mu$ g in a volume of  $\sim 3$   $\mu$ l) to be electroporated into sterile 1.5 ml microfuge tubes.
2. Thaw the competent cells at room temperature for several minutes. Add 50  $\mu$ l of cells to each DNA sample; gently pipet up and down to mix.
3. Incubate the samples at room temperature for 30 min.
4. Set the MicroPulser Electroporator to “StA.” See Section 4 for operating instructions.
5. Transfer the mixture of cells and DNA to a 0.2 cm electroporation cuvette and tap the suspension to the bottom of the tube. Place the cuvette in the chamber slide. Push the slide into the chamber until the cuvette is seated between the contacts in the base of the chamber.
6. Pulse once.
7. Remove the cuvette from the chamber and immediately add 1 ml of SMMP medium containing a subinhibitory concentration of antibiotic; gently transfer the cells to a sterile 17 x 100 mm tube using a Pasteur pipet. Incubate for 1 hr at 37°C, shaking at 250 rpm.
8. Check and record the pulse parameters. The time constant should be close to 2.5 milliseconds. The field strength can be calculated as actual kilovolts (kV) / cuvette gap (cm).
9. Plate aliquots of the electroporated cells on trypticase soy agar containing selective antibiotic. Incubate plates for 36–48 hr at 37°C.

## 6.3 Solutions and Reagents for Electroporation

1. B2 broth: 10 g casein hydrolysate, 25 g yeast extract, 5 g glucose, 25 g NaCl, 1 g  $\text{K}_2\text{HPO}_4$ ; dissolve in 900 ml water and adjust pH to 7.5; bring volume to 1.0 L. Autoclave.
2. SMMP: 55 ml 2x SMM, 40 ml 4x Penassay broth, 5 ml 10% (w/v) bovine albumin; adjust pH to 7.0; filter sterilize.
3. Trypticase soy agar: 40 g trypticase soy agar in 1.0 L of water. Autoclave.

4. 2x SMM: 25 ml 0.2 M sodium hydrogen maleate, 40 ml 0.1 N NaOH; adjust the pH to 6.5. Add 5 ml of 1 M MgCl<sub>2</sub>, 42.7 g sucrose; dissolve and bring volume to 125 ml. Filter sterilize.
5. 4x Penassay broth: 17.5 g Antibiotic Medium 3 dissolved in 250 ml water. Autoclave.
6. 0.2 M sodium hydrogen maleate: 11.6 g maleic anhydride or 13.7 g maleic acid, 4 g NaOH; dissolve in 500 ml water. Autoclave.

## Section 7

# Electroporation of *Agrobacterium tumefaciens*

## 7.1 Preparation of Electrocompetent Cells

See Lin 1995 for additional information.

1. Inoculate 1.5 L of YM broth in a 2.8 L Fernbach flask with an aliquot from log phase culture of *A. tumefaciens*.
2. Incubate at 30°C overnight, shaking at 300 rpm to a density of 5–10 × 10<sup>7</sup> cells/ml.
3. Decant the cells into sterile 500 ml centrifuge bottles and pellet the cells by centrifugation at 3,000 × g for 10 min at 4°C.
4. Carefully pour off and discard the supernatant; place the centrifuge bottles with the cell pellets on ice.
5. Add ~50 ml of sterile, ice-cold 10% glycerol to each of the bottles and vortex to resuspend the cell pellets; bring the volume in each of the centrifuge bottles to 500 ml with sterile, ice-cold 10% glycerol. Pellet the cells by centrifugation at 3,000 × g for 10 min at 4°C; pour off and discard the supernatant.
6. Wash the cells again as in step 5.
7. Resuspend each of the cell pellets in 5 ml of sterile, ice-cold 10% glycerol and transfer to a chilled 30 ml Oakridge tube. Pellet the cells by centrifugation at 3,000 × g for 5 min at 4°C; pour off and discard the supernatant.
8. Resuspend the cell pellet in 0.5 ml of sterile, ice-cold 1 M sorbitol; the final cell volume should be ~1.5 ml and the cell concentration should be ~5 × 10<sup>10</sup> cells/ml. Dispense 200 µl aliquots of the electrocompetent cells in sterile 1.5 ml microfuge tubes; freeze the cells in an isopropanol-dry ice bath, then store at –70°C. The cells are stable for about six months under these conditions.

## 7.2 Electroporation

1. Pipet the DNA samples (up to 5 µl) to be electroporated into sterile 1.5 ml microfuge tubes; the DNA should be in either water or TE. Place tubes on ice.
2. For each DNA sample to be electroporated, add 1 ml of YM broth to a 17 × 100 tube at room temperature, and place a 0.1 cm electroporation cuvette on ice.
3. Thaw the electrocompetent *A. tumefaciens* cells on ice. Add 20 µl of electrocompetent cells to each DNA sample to be electroporated; gently tap the tubes to mix.
4. Set the MicroPulser Electroporator to “Agr.” See Section 4 for operating instructions.

5. Transfer the DNA-cell samples to the electroporation cuvettes and tap the suspension to the bottom of the tube. Place the cuvette in the chamber slide. Push the slide into the chamber until the cuvette is seated between the contacts in the base of the chamber. Pulse once.
6. Remove the cuvette from the chamber and immediately use the YM broth in the 17 x 100 mm tube to transfer the cells from the cuvette to the tube.
7. Check and record the pulse parameters. The time constant should be about 5 milliseconds. The field strength can be calculated as actual kilovolts (kV) / cuvette gap (cm).
8. Incubate the cells for 3 hr at 30°C, shaking at 250 rpm. Plate aliquots of the electroporated cells on YM agar plates containing the appropriate selective medium. Incubate plates for 48 hr at 30°C.

### 7.3 Solutions and Reagents for Electroporation

1. YM broth: 0.4 g yeast extract, 10 g mannitol, 0.1 g NaCl, 0.1 g MgSO<sub>4</sub>, 0.5 g K<sub>2</sub>HPO<sub>4</sub>·3H<sub>2</sub>O; dissolve in 1.0 L of water and adjust pH to 7.0. Autoclave. For YM plates, add 15 g agar/1 L of YM broth.

## Section 8

# Electroporation of *Saccharomyces cerevisiae*

### 8.1 Preparation of Electrocompetent Cells

See Becker and Guarantee 1991 and Ausubel et al. 1987 for additional information.

1. Inoculate 500 ml of YPD in a 2.8 L Fernbach flask with an aliquot from an overnight culture of *S. cerevisiae*. The doubling time of *S. cerevisiae* is approximately 2 hr at 30°C.
2. Incubate at 30°C overnight, shaking at 250 rpm, to a density of ~1 x 10<sup>8</sup> cells/ml.
3. Chill the cells in an ice-water bath for 15 min to stop growth.
4. Decant the cells into two sterile 250 ml centrifuge bottles and pellet the cells by centrifugation at 3,000 x g for 5 min at 4°C.
5. Carefully pour off and discard the supernatant; place the centrifuge bottles with the cell pellets on ice.
6. Add ~50 ml of sterile, ice-cold water to each of the bottles and vortex to resuspend the cell pellets; bring the volume in each of the centrifuge bottles to 250 ml. Pellet the cells by centrifugation at 3,000 x g for 5 min at 4°C; pour off and discard the supernatant.
7. Wash the cells again as in step 6 with a total of 250 ml sterile, ice-cold water.
8. Resuspend the cell pellet in 20 ml of sterile, ice-cold 1 M sorbitol and transfer to a chilled 30 ml Oakridge tube. Pellet the cells by centrifugation at 3,000 x g for 5 min at 4°C; pour off and discard the supernatant.
9. Resuspend the cell pellet in 0.5 ml of sterile, ice-cold 1 M sorbitol; the final cell volume should be ~1.3 ml and the cell concentration should be ~1 x 10<sup>10</sup> cells/ml. Keep the cells on ice and use as soon as possible for electroporation.

## 8.2 Electroporation

1. Pipet the DNA samples (5–100 ng in a volume of 5  $\mu$ l) to be electroporated into sterile 1.5 ml microfuge tubes. Place tubes on ice.
2. If 0.2 cm cuvettes are used, add 40  $\mu$ l of the competent cells to each DNA sample; if 0.4 cm cuvettes are used, add 80  $\mu$ l of the competent cells to each DNA sample. Mix gently and incubate on ice for ~5 min.
3. Set the MicroPulser Electroporator to “Sc2” when using 0.2 cm cuvettes or to “Sc4” when using 0.4 cm cuvettes. See Section 4 for operating instructions.
4. Transfer the DNA-cell samples to the appropriate electroporation cuvettes that have been chilled in ice and tap the suspension to the bottom of the tube. Place the cuvette in the chamber slide. Push the slide into the chamber until the cuvette is seated between the contacts in the base of the chamber. Pulse once.
5. Remove the cuvette from the chamber and immediately add 1 ml of ice-cold 1 M sorbitol to the cuvette; gently transfer the diluted cells into a sterile tube.
6. Check and record the pulse parameters. The time constant should be close to 5 milliseconds. The field strength can be calculated as actual kilovolts (kV) / cuvette gap (cm).
7. Plate aliquots of the electroporated cells on selective agar plates containing 1 M sorbitol. Incubate plates for 48–72 hr at 30°C.

## 8.3 Solutions and Reagents for Electroporation

1. YPD: 10 g yeast extract, 20 g peptone, dissolve in 900 ml water. Autoclave. Add 100 ml sterile 20% glucose.
2. 1 M sorbitol: 182.2 g sorbitol, dissolve in 800 ml water. Bring volume to 1.0 L with water. Autoclave.
3. 20% glucose: 20 g glucose, dissolve in 60 ml water. Adjust volume to 100 ml with water. Sterilize through a 0.22  $\mu$ m filter.

# Section 9

## Electroporation of *Schizosaccharomyces pombe*

### 9.1 Preparation of Electrocompetent Cells

See Prentice 1991 for additional information.

1. Inoculate 500 ml of YCD in a 2.8 L Fernbach flask with an aliquot from an overnight culture of *S. pombe*. The doubling time of *S. pombe* is approximately 2 hr at 30°C.
2. Incubate at 30°C overnight, shaking at 250 rpm, to a density of  $1 \times 10^7$  cells/ml (OD 600 ~0.7).
3. Chill the cells in an ice-water bath for 15 min to stop growth.
4. Decant the cells into two sterile 250 ml centrifuge bottles and pellet the cells by centrifugation at 3,000 x g for 5 min at 4°C.

5. Carefully pour off and discard the supernatant; place the centrifuge bottles with the cell pellets on ice.
6. Add ~50 ml of sterile, ice-cold 1.2 M sorbitol to each of the bottles and vortex to resuspend the cell pellets; bring the volume in each of the centrifuge bottles to 250 ml. Pellet the cells by centrifugation at 3,000 x g for 5 min at 4°C; pour off and discard the supernatant.
7. Wash the cells again as in step 6 with a total of 250 ml sterile, ice-cold 1.2 M sorbitol.
8. Resuspend the cell pellet in 20 ml of sterile, ice-cold 1.2 M sorbitol and transfer to a chilled 30 ml Oakridge tube. Pellet the cells by centrifugation at 3,000 x g for 5 min at 4°C; pour off and discard the supernatant.
9. Resuspend the cell pellet in 0.5 ml of sterile, ice-cold 1.2 M sorbitol; the final cell volume should be ~1.3 ml and the cell concentration should be  $\sim 1 \times 10^9$  cells/ml. Keep the cells on ice and use as soon as possible for electroporation.

## 9.2 Electroporation

1. Pipet the DNA samples (up to 1  $\mu$ g) to be electroporated into sterile 1.5 ml microfuge tubes. Place tubes on ice.
2. Add 200  $\mu$ l of the competent cells to each DNA sample and mix gently.
3. Set the MicroPulser Electroporator to “ScS.” See Section 4 for operating instructions.
4. Transfer the DNA-cell samples to 0.2 cm electroporation cuvettes that have been chilled in ice and tap the suspension to the bottom of the tube. Place the cuvette in the chamber slide. Push the slide into the chamber until the cuvette is seated between the contacts in the base of the chamber. Pulse once.
5. Remove the cuvette from the chamber and immediately add 0.8 ml of ice-cold 1.2 M sorbitol to the cuvette; gently transfer the diluted cells to a sterile tube.
6. Check and record the pulse parameters. The time constant should be close to 5 milliseconds. The field strength can be calculated as actual kilovolts (kV) / cuvette gap (cm).
7. Incubate the tubes at room temperature for 40–60 min. Plate aliquots of the electroporated cells on minimal agar plates containing 1.2 M sorbitol. Incubate plates for 60–96 hr at 30°C.

## 9.3 Solutions and Reagents for Electroporation

1. YCD medium: 10 g yeast extract, 2 g casamino acids, dissolve in 900 ml water. Autoclave. Add 100 ml 20% glucose.
2. 1.2 M sorbitol: 218.6 g sorbitol, dissolve in 700 ml water. Add water to 1.0 L.

## Section 10

# Electroporation of *Dictyostelium discoideum*

### 10.1 Preparation of Electrocompetent Cells

See Howard et al. 1988 and Knecht and Pang 1995 for additional information.

1. Inoculate *D. discoideum* cells at a concentration of  $5\text{--}7 \times 10^5$  cells/ml into 40 ml of HL5 medium in a 500 ml flask. The cells may either be scraped from a plate or transferred from liquid medium. The doubling time of *D. discoideum* is ~12 hr at 21°C.
2. Incubate the culture at 21°C for about 24 hr, shaking at 125 rpm. About 16–20 hr prior to preparing the competent cells, dilute the cells to  $7 \times 10^5$  cells/ml with HL5 medium. Incubate at 21°C overnight, shaking at 125 rpm.
3. Transfer 100 ml of the cells into two sterile, disposable, 50 ml centrifuge tubes and incubate on ice for 15 min to stop growth.
4. Pellet the cells by centrifugation at 400 x g for 5–7 min at room temperature.
5. Carefully pour off and discard the supernatant; place the centrifuge bottles with the cell pellets on ice.
6. Pool the cell pellets and resuspend in 50 ml of sterile, ice-cold E buffer. Pellet the cells by centrifugation at 400 x g for 5–7 min at room temperature.
7. Carefully pour off and discard the supernatant; place the centrifuge bottles with the cell pellets on ice and resuspend the cells at a concentration of  $1 \times 10^7$  cells/ml. Keep the cells on ice and use as soon as possible for electroporation.

### 10.2 Electroporation

1. Pipet the DNA samples (up to 50 µg) to be electroporated into sterile 1.5 ml microfuge tubes. Place tubes on ice.
2. Add 800 µl of the competent cells to each DNA sample and pipet up and down to mix; incubate on ice for ~1 min.
3. Set the MicroPulser Electroporator to “Dic.” See Section 4 for operating instructions.
4. Transfer the DNA-cell samples to 0.4 cm electroporation cuvettes that have been chilled on ice and tap the suspension to the bottom of the tube. Place the cuvette in the chamber slide. Push the slide into the chamber until the cuvette is seated between the contacts in the base of the chamber. Pulse once (the program delivers two pulses approximately 5 seconds apart).
5. Remove the cuvette from the chamber and immediately dilute the cells to 10 ml with the appropriate medium.
6. Check and record the pulse parameters. The time constant should be 1 millisecond. The field strength can be calculated as actual kilovolts (kV) / cuvette gap (cm).
7. When selecting for complementation of an auxotrophic mutant, the cells may be plated immediately into selective medium lacking the appropriate nutrient. When selecting for antibiotic resistance, incubate the cells overnight at 21°C prior to adding the selective agent.

### 10.3 Solutions and Reagents for Electroporation

1. HL5 medium: 17.8 g bacteriological peptone (Oxoid, Ogdensburg, NY), 7.2 g yeast extract, 0.54 g Na<sub>2</sub>HPO<sub>4</sub>, 0.4 g KH<sub>2</sub>PO<sub>4</sub>, 130 µl B12/folic acid mix; bring to 1.0 L with water and adjust pH to 6.3–6.5. Autoclave for 25 min on two successive days. Prior to use, add 20 ml of 50% glucose and 10 ml of 100 x Antibiotic-Antimycotic (Thermo Fisher Scientific Inc.).
2. B12/folic acid mix: 5 mg B12, 200 mg folic acid; add 95 ml water, then adjust pH to 6.5–6.8 with 5N NaOH; bring to 100 ml with water. Filter sterilize and store at –20°C protected from light.
3. E buffer: 10 ml 100 mM NaH<sub>2</sub>PO<sub>4</sub> adjusted to pH 6.1 with KOH, 10 ml 0.5 M sucrose, 80 ml water; autoclave.

## Section 11

### Electroporation of *Pichia pastoris*

#### 11.1 Preparation of Electrocompetent Cells

See Cregg and Russell 1998 for additional information.

1. Inoculate 500 ml of YPD in a 2.8 L Fernbach flask with an aliquot from a fresh overnight culture of *P. pastoris*. The doubling time of *P. pastoris* is approximately 2 hr at 30°C.
2. Incubate at 30°C overnight, shaking at 300 rpm, to a density of 5–7 x 10<sup>7</sup> cells/ml.
3. Decant the cells into two sterile 250 ml centrifuge bottles and pellet the cells by centrifugation at 3,000 x g for 5 min at 4°C.
4. Carefully pour off and discard the supernatant.
5. Add 50 ml of sterile YPD/HEPES to each of the bottles and vortex to resuspend the cell pellets; add 1.25 ml of 1 M DTT to each bottle; mix gently. Incubate the cells for 15 min at 30°C.
6. Add 200 ml of sterile, ice-cold 1 M sorbitol to each centrifuge bottle. Pellet the cells by centrifugation at 3,000 x g for 5 min at 4°C; pour off and discard the supernatant.
7. Add ~50 ml of sterile, ice-cold 1 M sorbitol to each of the bottles and vortex to resuspend the cell pellets; bring the volume in each of the centrifuge bottles to 250 ml with sterile, ice-cold 1 M sorbitol. Pellet the cells by centrifugation at 3,000 x g for 5 min at 4°C; pour off and discard the supernatant.
8. Resuspend each cell pellet in 10 ml of sterile, ice-cold 1 M sorbitol and pool in a chilled 30 ml Oakridge tube. Pellet the cells by centrifugation at 3,000 x g for 5 min at 4°C; pour off and discard the supernatant.
9. Resuspend the cell pellet in 0.5 ml of sterile, ice-cold 1 M sorbitol; the final cell volume should be ~1.3 ml and the cell concentration should be ~1 x 10<sup>9</sup> cells/ml. Keep the cells on ice and use as soon as possible for electroporation.

## 11.2 Electroporation

1. Pipet the DNA samples (up to 10 µg) to be electroporated into sterile 1.5 ml microfuge tubes. Place tubes on ice.
2. Add 40 µl of the competent cells to each DNA sample and mix gently.
3. Set the MicroPulser Electroporator to "Pic." See Section 4 for operating instructions.
4. Transfer the DNA-cell samples to 0.2 cm electroporation cuvettes that have been chilled in ice and tap the suspension to the bottom of the tube. Place the cuvette in the chamber slide. Push the slide into the chamber until the cuvette is seated between the contacts in the base of the chamber. Pulse once.
5. Remove the cuvette from the chamber and immediately add 1.0 ml of ice-cold 1 M sorbitol to the cuvette when selecting for complementation of an auxotrophic mutant or 1.0 ml of ice-cold YPD/sorbitol when selecting for antibiotic resistance. Gently transfer the diluted cells to a sterile tube.
6. Check and record the pulse parameters. The time constant should be close to 5 milliseconds. The field strength can be calculated as actual kilovolts (kV) / cuvette gap (cm).
7. When selecting for complementation of an auxotrophic mutant, the cells may be plated immediately onto minimal agar plates containing 1 M sorbitol but lacking the appropriate nutrient. When selecting for antibiotic resistance, incubate the cells at 30°C for 1–2 hr without shaking; plate aliquots of the electroporated cells on YPD agar plates containing 1 M sorbitol with the appropriate antibiotic. Incubate the plates for 72–96 hr at 30°C.

## 11.3 Solutions and Reagents for Electroporation

1. YPD/HEPES: 100 ml YPD medium, 20 ml 1 M HEPES, pH 8.0.
2. 1 M DTT: 1.55 g dithiothreitol, dissolve in 8 ml water. Bring the volume to 10 ml with water. Filter sterilize.
3. YPD/sorbitol: 10 g yeast extract, 20 g peptone, 182.2 g sorbitol, dissolve in 700 ml water; bring volume to 900 ml with water. Autoclave. Add 100 ml sterile 20% glucose.



# Appendix I

## References

- Allen SP and Blaschek HP (1990). Factors involved in the electroporation-induced transformation of *Clostridium perfringens*. FEMS Microbiol Lett 70, 217–220.
- Ausubel FM et al. (1987). Current Protocols in Molecular Biology (New York: John Wiley & Sons).
- Becker DM and Guarente L (1991). High-efficiency transformation of yeast by electroporation. Methods Enzymol 194, 182–187.
- Böttger EC (1988). High-efficiency generation of plasmid cDNA libraries using electro-transformation. Biotechniques 6, 878–880.
- Chang DC et al. (1991). Guide to Electroporation and Electrofusion (San Diego: Academic Press Inc.).
- Cregg JM and Russell KA (1998). Transformation. Methods Mol Biol 103, 27–39.
- Cruz-Rodz AL and Gilmore MS (1990). High efficiency introduction of plasmid DNA into glycine treated *Enterococcus faecalis* by electroporation. Mol Gen Genet 224, 152–154.
- Dennis JJ and Sokol PA (1995). Electrotransformation of *Pseudomonas*. Methods Mol Biol 47, 125–133.
- Dower WJ (1990). Electroporation of bacteria: A general approach to genetic transformation. In Genetic Engineering — Principles and Methods, J. K. Setlow, ed. (New York: Plenum Publishing Corp.), pp. 275–295.
- Dower WJ et al. (1988). High efficiency transformation of *E. coli* by high voltage electroporation. Nucleic Acids Res 16, 6,127–6,145.
- Dower WJ et al. (1992). Protocols for the transformation of bacteria by electroporation. In Guide to Electroporation and Electrofusion, D. C. Change et al., eds. (San Diego: Academic Press Inc.), p. 485.
- Grey M and Brendel MF (1995). Ten-minute electrotransformation of *Saccharomyces cerevisiae*. In Methods in Molecular Biology, J. A. Nickoloff, ed. (Totowa, NJ: Humana Press) 47, pp. 269–272.
- Heery DM and Dunican LK (1989). Improved efficiency M13 cloning using electroporation. Nucleic Acids Res 17, 8,006.
- Holo H and Nes (1995). IF Transformation of *Lactococcus* by electroporation. Methods Mol Biol 47, 195–199.
- Howard PK et al. (1988). Establishment of a transient expression system for *Dictyostelium discoideum*. Nucleic Acids Res 16, 2,613–2,623.
- Jacobs M et al. (1990). High-efficiency electro-transformation of *Escherichia coli* with DNA from ligation mixtures. Nucleic Acids Res 18, 1,653.
- Knecht D and Pang KM (1995). Electroporation of *Dictyostelium discoideum*. Methods Mol Biol 47, 321–330.
- Lee JC (1995). Electrotransformation of staphylococci. Methods Mol Biol 47, 209–216.
- Leonardo ED and Sedivy JM (1990). A new vector for cloning large eukaryotic DNA segments in *Escherichia coli*. Biotechnology 8, 841–844.
- Lin J-J (1995). Electrotransformation of *Agrobacterium*. Methods Mol Biol 47, 171–178.
- Miller EM and Nickoloff JA (1995). *Escherichia coli* electrotransformation. Methods Mol Biol 47, 105–113.
- Park SF and Stewart GS (1990). High efficiency transformation of *Listeria monocytogenes* by electroporation of penicillin-treated cells. Gene 94, 129–132.
- Prentice HL (1992). High efficiency transformation of *Schizosaccharomyces pombe* by electroporation. Nucleic Acids Res 20, 621.
- Shigekawa K and Dower WJ (1988). Electroporation of eukaryotes and prokaryotes: A general approach to the introduction of macromolecules into cells. Biotechniques 6, 742–751.
- Siguret V et al. (1994). Effect of plasmid size on transformation efficiency by electroporation of *Escherichia coli* DH5 alpha. Biotechniques 16, 422–426.
- Simon D and Ferretti JJ (1991). Electrotransformation of *Streptococcus pyogenes* with plasmid and linear DNA. FEMS Microbiol Lett 66, 219–224.
- Somkuti GA and Steinberg DH (1988). Genetic transformation of *Streptococcus thermophilus* by electroporation. Biochimie 70, 579–585.
- Summers DK and Withers HL (1990). Electrotransfer: Direct transfer of bacterial plasmid DNA by electroporation. Nucleic Acids Res 18, 2192.
- Taketo A (1988). DNA transfection of *Escherichia coli* by electroporation. Biochim Biophys Acta 949, 318–324.
- Taketo A (1989). RNA transfection of *Escherichia coli* by electroporation. Biochim Biophys Acta 1,007, 127–129.
- Trevors JT (1992). Electrotransformation of bacteria by plasmid DNA. In Guide to Electroporation and Electrofusion, Chang D. C. et al. eds. (San Diego: Academic Press Inc.), p. 265.
- Willson TA and Gough NM (1988). High voltage *E. coli* electro-transformation with DNA following ligation. Nucleic Acids Res 16, 11,820.
- Zabarovsky ER and Winberg G (1990). High efficiency electroporation of ligated DNA into bacteria. Nucleic Acids Res 18, 5,912.

## Appendix II

# Troubleshooting Guide for the MicroPulser Electroporator

Operational problem	Possible cause and solution
1. Display does not light when unit is turned on.	<ul style="list-style-type: none"> <li>■ Power is not supplied to electronics. Check power cord and wall outlet power source</li> <li>■ Check that power switch is on</li> <li>■ Check/replace fuse. Two 2 A, 250 V Type T fuses are located on the back of the instrument immediately above the power cord</li> </ul>
2. When the buttons are pressed, the unit does not indicate “PLS.”	<ul style="list-style-type: none"> <li>■ No pulse delivery</li> <li>■ Pulse button is not depressed hard enough</li> <li>■ Verify instrument operation by the following tests:               <ul style="list-style-type: none"> <li>– Turn on the power switch</li> <li>– Verify the LED next to “Bacteria” is illuminated. Press the “Settings” button</li> <li>– Verify the LED next to “Fungi” is illuminated. Press the “Settings” button again</li> <li>– Verify the LED next to “Manual” is illuminated. Press the “Settings” button again</li> <li>– The LED next to “Bacteria” should be lit. Press the “Raise” button several times to change the display from “Ec1” to “Ec3.” Press the “Raise” and “Lower” buttons simultaneously to verify the display reads “3.00” volts</li> <li>– Press the “Pulse” button until “PLS” is displayed; a beep tone should be heard within 6 sec</li> <li>– Press the “Measurements” button. Verify the “Actual kV” LED illuminates. Verify the display LED shows a reading between “2.95” and “3.05”</li> <li>– Press the “Measurements” button again. Verify the “Time ms” LED illuminates</li> <li>– Verify the display LED shows a reading between “5.50” and “6.50.” If problems persist, contact Bio-Rad</li> </ul> </li> </ul>

Electrical problem	Possible cause and solution
1. Instrument shows “Arc” on display LED.	<p data-bbox="621 243 1417 478">Arcing in the cuvette is the result either of an actual arc occurring or of medium that is too conductive. An actual arc will occur usually only at high voltage (&gt;1,500 V). If a medium is too conductive, the display LED may show “Arc” even though an actual arc has not occurred. The limit of conductivity depends on the voltage, electrode gap, and sample volume, but under standard conditions, solutions of 10 meq or higher (&lt;600 <math>\Omega</math> resistance) will certainly indicate “Arc.”</p> <p data-bbox="621 510 1417 745">To determine whether an actual arc has occurred or the medium is too conductive, look at the cuvette while pulsing the sample again. If a small spark occurs across the electrodes, an arc has occurred. If no spark is observed, lower the voltage and repulse the sample. Continue lowering and repulsing the sample until a pulse occurs. Check the time constant; if the time constant is low (&lt;3 ms), the sample is too conductive.</p> <p data-bbox="621 756 1417 793">There are several causes of excessive conductivity:</p> <ol data-bbox="621 804 1417 1182" style="list-style-type: none"> <li data-bbox="621 804 1417 877">1. Washing and resuspending cells in a buffer too high in ionic strength.</li> <li data-bbox="621 888 1417 982">2. Insufficient washing of the cells — salts from the growth medium are not completely removed; the cells should be washed at least three to four times with nonconductive solution.</li> <li data-bbox="621 993 1417 1066">3. Lysed cells in the preparation — cell contents contribute to conductivity.</li> <li data-bbox="621 1077 1417 1182">4. DNA solution too high in salt concentration; for example, CsCl carried over from plasmid preparation, or residual salts from ethanol precipitation or ligation.</li> </ol> <p data-bbox="621 1192 1417 1266">Electroporation with cuvettes above 0°C: reducing sample temperature increases sample resistance.</p>
2. Wrong time constant.	<p data-bbox="621 1276 1417 1482">Samples electroporated in the MicroPulser Electroporator should have a time constant close to 5 ms. If the time constant is much shorter than the expected value (e.g., 3 ms instead of 5 ms), the sample is too conductive. The probable reasons for this are listed above under “arcing.” Correct the problem of high conductivity by additional washing of the cells, or by removal of salts from the DNA preparation.</p>
3. Sample does not “twitch.”	<p data-bbox="621 1493 1417 1629">This may mean that the pulse is not reaching the sample. Check the connections between the MicroPulser Electroporator and sample chamber. Check to see that the contacts in the base of the sample chamber are not broken.</p>
4. Instrument displays “no” on front panel.	<p data-bbox="621 1640 1417 1707">The manual setting is used and the voltage is set to 0.00. Use the “Raise” button to select a voltage between 0.02 and 3.00 kV.</p>
5. Instrument displays “err” on front panel.	<p data-bbox="621 1717 1417 1780">Turn instrument off, then on again. If problem persists, contact Bio-Rad.</p>

---

## Biological

---

The general symptom addressed in this section is transformation efficiencies that are too low to detect or too low to be useful. The following is a list of the areas of possible problems and some suggested solutions.

Problem	Possible cause and solution
1. The pulse.	<p>Is the pulse applied to the sample? At high voltage with a small-volume (40 <math>\mu</math>l) sample this is easy to check. The sample will “twitch” when pulsed. If you don’t see a twitch, refer to the electrical troubleshooting section for information on electrical problems. Also, make sure that the cuvette is in contact with the electrodes at the back of the sample chamber. If electrodes are broken or corroded, call Bio-Rad for replacements.</p> <p>Are the amplitude and length of the pulse sufficient? <i>E. coli</i> requires pulses of approximately 5 ms with field strengths of 12 to 18 kV/cm. <i>S. cerevisiae</i> requires pulses of approximately 5 ms with a field strength of ~7.5 kV/cm. There is usually some cell death with electrical conditions producing transformation. Survival rates of 20 to 80% are typical. If no cell death occurs, the pulse is probably too weak. Conversely, if too many cells are killed (&gt;80%), the pulse is probably too intense and transformation will probably be poor. To find the optimum pulse characteristics, use a pulse length of ~5 ms and test for transformation over a range of field strengths.</p>
2. The DNA.	<p>Check the quantity and quality of the DNA on a gel. Often, minipreps contain less DNA than expected. DNA stored improperly for long periods may be degraded and may lack transforming activity.</p> <p>Some preparations of DNA may contain substances that inhibit transformation or are toxic to the cells. Try to use DNA free of SDS, phenol, etc.</p> <p>Is the selection appropriate for the marker (and its level of expression)?</p>
3. The cells.	<p>Were the cells harvested at the correct stage in the growth phase? Bacterial cells generally show the highest transformation efficiencies when harvested in the early- to mid-log growth phase. Yeast cells generally show the highest transformation efficiencies when harvested in late-log phase. Different growth conditions may improve transformation.</p> <p>Are too many cells killed? The pulse is too intense, toxic substances are present in DNA or cell preparations, or wrong temperature of electroporation are all possibilities.</p> <p>Are the cells transferred to outgrowth medium immediately after the pulse? For <i>E. coli</i> this is very important.</p> <p>Is the correct selection applied after the recovery period?</p>
4. The temperature.	<p>Are the cuvettes cold?</p> <p>Is the cuvette holder (slide) prechilled?</p> <p>If frozen, have the cells been stored properly (usually in 10–15% glycerol at <math>-70^{\circ}\text{C}</math>)?</p>

---

## Appendix III

# Product Information

### Specifications

Input voltage	100–120 V RMS, 50/60 Hz; automatic mains voltage switching 220–240 V RMS, 50/60 Hz
Input current	2 A RMS (100–120 V), 1 A RMS (220–240 V)
Maximum output voltage and current	3,000 V peak into >3.3 k $\Omega$ load limited to 100 A peak maximum
Output waveform	Decaying exponential waveform with RC time constant of 5 ms, assuming loads of ~3.3 k $\Omega$
Output voltage adjustment	Voltage adjustable in 200–3,000 V range with 10 V display resolution; 10 preprogrammed voltage steps
Dimensions (H x W x D)	8 x 21 x 31 cm
Weight	2.9 kg (6.4 lb)

### Environmental Conditions

Operating temperature	3.5–35°C, indoor use
Humidity	0–95%, in the absence of condensation
Altitude	Up to 2,000 m
Pollution degree	2

### Input Power

Supply voltage fluctuations	$\pm 10\%$ (unless otherwise specified) Maximum current 2 A Maximum power 240 W
Overvoltage category	II

## Related Products

### Cuvettes

Catalog # Description

#### Mini Packs

- 1652083 **Gene Pulser/MicroPulser Electroporation Cuvettes**, 0.1 cm gap, 5 sterile
- 1652082 **Gene Pulser/MicroPulser Electroporation Cuvettes**, 0.2 cm gap, 5 sterile
- 1652081 **Gene Pulser/MicroPulser Electroporation Cuvettes**, 0.4 cm gap, 5 sterile

#### Standard Packs

- 1652089 **Gene Pulser/MicroPulser Electroporation Cuvettes**, 0.1 cm gap, 50 sterile
- 1652086 **Gene Pulser/MicroPulser Electroporation Cuvettes**, 0.2 cm gap, 50 sterile
- 1652088 **Gene Pulser/MicroPulser Electroporation Cuvettes**, 0.4 cm gap, 50 sterile

#### Jumbo Packs

- 1652093 **Gene Pulser/MicroPulser Electroporation Cuvettes**, 0.1 cm gap, 500 sterile
- 1652092 **Gene Pulser/MicroPulser Electroporation Cuvettes**, 0.2 cm gap, 500 sterile
- 1652091 **Gene Pulser/MicroPulser Electroporation Cuvettes**, 0.4 cm gap, 500 sterile

### Plasmid Purification

Catalog # Description

- 7326100 **Quantum Prep Plasmid Miniprep Kit**, 100 preps, plasmid DNA purification kit, includes Quantum Prep matrix, solutions (cell resuspension and lysis, neutralization, wash), spin columns
- 7326120 **Quantum Prep Plasmid Midiprep Kit**, 20 preps, large quantity plasmid DNA purification kit, includes Quantum Prep matrix, solutions (cell resuspension and lysis, neutralization, wash), midi spin columns



Visit [bio-rad.com/MicroPulsar](http://bio-rad.com/MicroPulsar) for more information.

BIO-RAD, GENE PULSER, and QUANTUM PREP are trademarks of Bio-Rad Laboratories, Inc. in certain jurisdictions. All trademarks used herein are the property of their respective owner.



**Bio-Rad  
Laboratories, Inc.**

Life Science  
Group

---

**Website** [bio-rad.com](http://bio-rad.com) **USA** 1 800 424 6723 **Australia** 61 2 9914 2800 **Austria** 00 800 00 24 67 23 **Belgium** 00 800 00 24 67 23 **Brazil** 4003 0399 **Canada** 1 905 364 3435 **China** 86 21 6169 8500 **Czech Republic** 00 800 00 24 67 23 **Denmark** 00 800 00 24 67 23 **Finland** 00 800 00 24 67 23 **France** 00 800 00 24 67 23 **Germany** 00 800 00 24 67 23 **Hong Kong** 852 2789 3300 **Hungary** 00 800 00 24 67 23 **India** 91 124 4029300 **Israel** 0 3 9636050 **Italy** 00 800 00 24 67 23 **Japan** 81 3 6361 7000 **Korea** 82 2 3473 4460 **Luxembourg** 00 800 00 24 67 23 **Mexico** 52 555 488 7670 **The Netherlands** 00 800 00 24 67 23 **New Zealand** 64 9 415 2280 **Norway** 00 800 00 24 67 23 **Poland** 00 800 00 24 67 23 **Portugal** 00 800 00 24 67 23 **Russian Federation** 00 800 00 24 67 23 **Singapore** 65 6415 3188 **South Africa** 00 800 00 24 67 23 **Spain** 00 800 00 24 67 23 **Sweden** 00 800 00 24 67 23 **Switzerland** 00 800 00 24 67 23 **Taiwan** 886 2 2578 7189 **Thailand** 66 2 651 8311 **United Arab Emirates** 36 1 459 6150 **United Kingdom** 00 800 00 24 67 23

---

