

Instructions to update CFX Opus (Dx and RUO) Instrument Software

Step 1: Copy to the root directory of a USB drive the zipped file: Setup_3G-brelease-v1.5.144.0505_BT-RC-01_gca8333cd_UpgradeE

IMPORTANT: Do not unzip the file.

Step 2: Insert the USB drive on the CFX Opus instrument.

Step 3: Power on the CFX Opus instrument and log in as ADMIN.

NOTE: Only a user with an Admin role can perform the software update.

Step 4: From the Home screen, navigate to the Tools menu and select System Update.

Step 5: While the upgrade is in progress, do not shut down the system, and do not remove the USB drive. The CFX Opus instrument will automatically restart when the update is complete.